

# Model MKFT 115 | Dynamic climate chambers for rapid temperature changes with humidity control and extended low temperature range

The BINDER MKFT series environmental simulation chamber for low temperature alternating climate profiles is the specialist for dynamic alternating climate change between -70 °C and 180 °C. Large performance reserves and rapid cooling make the MKFT environmental simulation chamber the high-end product for complex standards-based climate tests.

# BENEFITS

- Homogeneous climate conditions thanks to APT.line™ technology
- Automatic water and wastewater management
- Pressure humidification with fast response times
- Comprehensive programming and data acquisition
- Large heated viewing window



Model 115

## **MAIN FEATURES**

- Temperature range: -70 °C to +180 °C
- Humidity range: 10 % to 98 % RH
- Integrated water-storage tank, 20 liters
- 4 zero-voltage relay contacts
- APT.line<sup>™</sup> preheating chamber technology
- Programmable condensation protection for test material
- Intuitive touchscreen controller with time-segment and real-time programming
- Internal data logger, measured values can be read out in open format via USB
- Heated viewing window with LED interior lighting
- Humidity regulation with capacitative humidity sensor and vapor humidification
- BINDER Multi Management Software APT-COM™ Basic Edition
- Troubleshooting system with visual and audible alarms
- Class 2 independent adjustable temperature safety device (DIN 12880) with visual alarm
- Access port with silicone plug: 50 mm, left



Model 115

- 4 castors, two with brakes
- Computer interface: Ethernet
- Adjustable ramp function
- Integrated chart recorder
- Real-time clock
- Door heating
- Alarm notification in the event of insufficient water in fresh water tank
- 1 stainless steel rack
- Complete safety connection kit for water supply and drainage, up to 1 m in height
- Inner chamber made of stainless steel
- CFC-free refrigerants R-452A and R-23
- Cooling with cascade compressor cooling unit

### **ORDERING INFORMATION**

Interior volume [L]	Voltage	Option model	Version	ArtNo.
	400 V 3~ ph 50 Hz	Standard	MKFT115-400V	9020-0382
115	480 V 3~ ph 60 Hz	with voltage and frequency converter	MKFT115-400V MKFT115-480V-C	9020-0362




# **TECHNICAL DATA**

Description	MKFT115-400V <sup>1</sup>	MKFT115-480V-C <sup>1</sup>
Article Number	9020-0382	9020-0362
Performance Data Temperature		
Temperature range	-70180	-70180
Temperature variation depending on setpoint [± K]	0.11.3	0.11.3
Temperature fluctuation depending on setpoint [± K]	0.10.1	0.10.1
Average heating-up rate according to IEC 60068-3-5 [K/min]	5.5	5.5
Average cooling down time according to IEC 60068-3-5 [K/min]	4.2	4.2
Max. heat compensation at 25 °C [W]	1500	1500
Performance Data Climate		
Temperature range [°C]	1095	1095
Temperature fluctuation depending on setpoint [± K]	0.11	0.11
Humidity range	1098	1098
Humidity fluctuation depending on setpoint	≤2,5 ± % RH	≤2,5 ± % RH
Dew point temperature range [°C]	594	594
Electrical data		
Rated Voltage [V]	400	480
Power frequency [Hz]	50	60
Nominal power [kW]	5	5
Unit fuse [A]	16	16
Phase (Nominal voltage)	3~	3~
Measures		
Interior volume [L]	115	115
Net weight of the unit (empty) [kg]	330	494
Permitted load [kg]	60	60
Load per rack [kg]	30	30
Viewing window width [mm]	288	288
Viewing window height [mm]	222	222
Wall clearance back [mm]	300	300
Wall clearance sidewise [mm]	200	200
Internal Dimensions		
Width [mm]	600	600
Height [mm]	480	480
Depth [mm]	400	400
Doors		
Unit doors	1	1
Housing dimensions not incl. fittings and connections		
Width net [mm]	980	980
Height net [mm]	1725	1725
Depth net [mm]	865	865
Environment-specific data		
Sound-pressure level [dB(A)]	64	67
Fixtures		
Number of shelves (std./max.)	1/4	1/4


1 All technical data is specified for unloaded units with standard equipment at an ambient temperature of +22 °C ±3 °C and a power supply voltage fluctuation of ±10 %. The temperature data is determined in accordance to BINDER factory standard following DIN 12880, observing the recommended wall clearances of 10 % of the height, width, and depth of the inner chamber. Technical data refers to 100 % fan speed. All indications are average values, typical for units produced in series. We reserve the right to change technical specifications at any time.

BINDER


# DIMENSIONS incl. fittings and connections [mm]

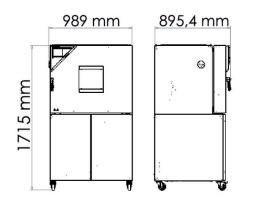


# CHARTS

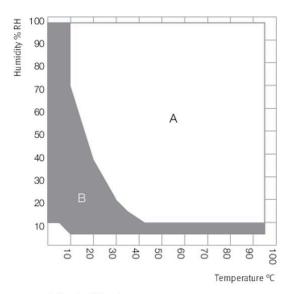


A: Standard Climate range B: Extended range with dry-air purge

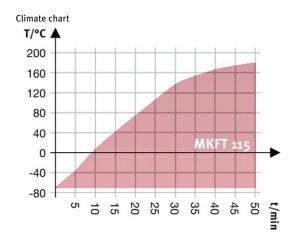



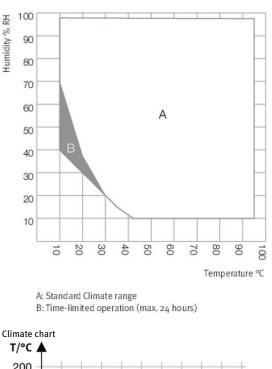


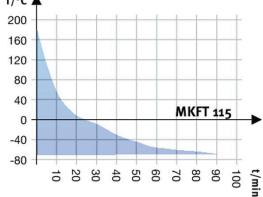

Cooling down rate

Humidity % RH



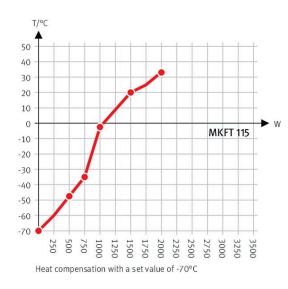

Heat compensation chart

#### **OPTIONS**

Description	*	ArtNo.
left		
30 mm	01	8012-0509
50 mm	01	8012-0511
80 mm	01	8012-0513
100 mm	01	8012-0836
125 mm	01	8012-0533
right		
30 mm	01	8012-0508
50 mm	01	8012-0510
8o mm	01	8012-0512
100 mm	01	8012-0835
125 mm	01	8012-0532
top		
8o mm	01	8012-0834
100 mm	01	8012-0832
125 mm	01	8012-0833
for temperature and humidity values (output not adjustable)	-	8012-1786
for temperature and humidity; for extending the measurement in center of chamber to include another test value	-	8012-1193
temperature measurement incl. certificate and 27 measuring points at specified temperature	-	8012-1608
temperature measurement incl. certificate, 15- 18 measuring points at specified temperature	-	8012-1588
temperature measurement incl. certificate, 9 measuring points at specified temperature	-	8012-1567
Measurement in center of chamber at 25 °C / 60% RH or at specified test values	-	8012-1187
temperature (according to DIN12880) and humidity measurement incl. certificate, 27 temperature measuring points and 1 humidity measuring point, at 25 °C / 60 % RH or at specified values	-	8012-1614
for the connection to an existing pressurized air network	-	8012-1803
lockable door handle	-	8012-1783
controlled, incl. connection; suitable for the simulation of current automobile industry standards	-	8012-1789
	left 30 mm 50 mm 80 mm 100 mm 125 mm right 30 mm 50 mm 80 mm 100 mm 125 mm 100 mm	Descriptionleft30 mm0130 mm0150 mm01100 mm01125 mm01right0130 mm0150 mm0180 mm01100 mm01100 mm01100 mm01125 mm01100 mm01125 mm01top0180 mm01100 mm01125 mm01for temperature and humidity values (output not adjustable)-for temperature and humidity; for extending the measurement in center of chamber to include another test value-temperature measurement incl. certificate and 27 measuring points at specified temperature-temperature measurement incl. certificate, 9 measuring points at specified temperature-Measurement in center of chamber at 25 °C / 60 % RH or at specified test values-temperature (according to DIN12880) and humidity measurement incl. certificate, 27 temperature measuring points and 1 thumidity measuring points at specified values-temperature (according to DIN12880) and humidity measurement incl. certificate, 27 temperature measuring points and 1 thumidity measuring points at specified values-temperature (according to DIN12880) and humidity measurement incl. certificate, 27 temperature measuring points and 1 thumidity measuring points at specified values-for the connection to an existing pressurized air network-

\* Notes > See last page



Designation	Description	*	ArtNo.
Pt 100 temperature sensor	additional flexible Pt 100, interior, for displaying the temperature on the unit display	-	8012-1791
RS 485 interface, 2-wire	Additional serial interface can be used parallel to Ethernet, for Multi Management Software APT-COM™	-	8012-1767
Temperature safety device class 2	with visual alarm (DIN 12880)	-	8012-1793
Water circuit	allows condensed water to be reused	-	8012-0785

\* Notes > See last page

### ACCESSORIES

Designation	Description	*	ArtNo.
APT-COM™ 4 GLP- Edition	for working under GLP-compliant conditions. Measured values are documented in a tamper-proof way in line with the requirements of FDA Regulation 21 CFR 11.		
	version 4, GLP edition	-	9053-0042
APT-COM™ 4	convenient unit and user management built on the BASIC edition. Suitable for networking up to 100 units.		
PROFESSIONAL-Edition	version 4, PROFESSIONAL edition	-	9053-0040
BINDER PURE AQUA SERVICE	System for preparation or complete desalination of tap water, complete set containing PURE AQUA 300 single-use cartridge, measuring device, and all necessary connecting parts	-	8012-0759
BINDER PURE AQUA SERVICE, accessories	Single-use, replacement cartridge for BINDER PURE AQUA System	-	6011-0165
pH-neutral detergent	concentrated, for gentle remove of residual contaminants; 1 kg	-	1002-0016
	IQ/OQ documents – supporting documents for validation performed by customers, consisting of: IQ/OQ checklists incl. calibration guide and comprehensive unit documentation; parameters: temperature and humidity values		
	Digital in PDF format	-	7057-0002
Qualification documents	Hard copy inside folder	-	7007-0002
Qualification documents	IQ/OQ/PQ documents – supporting documents for validation performed by customers, according to customer requirements, PQ section added to qualification folder IQ/OQ; parameters: temperature and humidity values		
	Digital in PDF format	-	7057-0006
	Hard copy inside folder	-	7007-0006
Rack	stainless steel	-	8012-2048
Rack accessories	fasteners (1 set of 4) for additional security of racks	-	8012-0620
Rack, reinforced	stainless steel, with fasteners (1 set of 4)	-	8012-0709
	RS 422 cable set and RS 485 / RS 422 interface converter for connection to 10-way plug distributor		
RS 485 / RS 422 interface converter	115 V option model	-	8012-0599
	230 V option model	-	8012-0589
Shelf, perforated	Stainless steel	-	6004-0030
WLAN kit	The kit contains a client bridge that establishes a wireless connection between BINDER units and APT-COM4, LIMS, or customer-specific software via an Ethernet interface. It provides an alternative solution in situations where units need to be positioned in locations without an on-site Ethernet connection. In secured networks, installation and configuration must be performed by the customer's IT service.	-	8500-0164

\* Notes > See last page

### SERVICES

Designation	Description	*	ArtNo.
Installation services			
Unit commissioning	Connect the unit to the customer-side connections (electricity, water, wastewater, gas), basic functions check, brief operating instructions. (excl.: unpacking, setup, controller instructions, programming, installation work)	05	DL10-0300
Unit instructions	Instruction regarding operating principle and basic functions of the unit, operation of the control electronics including programming	05	DL10-0700
Maintenance contracts			
Maintenance services			
BRONZE 3-year maintenance contract	Maintenance service as contractually agreed, visual inspection of mechanical and electrical components, check of control response, 20% discount on spare parts	05	DL20-0710
* Notes › See last page			

Data Sheet Model MKFT 115

# BINDER

Designation	Description	*	ArtNo.
GOLD 3-year maintenance contract	Maintenance service as contractually agreed, visual inspection of mechanical and electrical components, check of control response, 20% discount on spare parts, testing of all key functions, replacement of wear parts, calibration of one temperature and CO <sub>2</sub> value, including certificate	05	DL20-0940
Leakage test	Testing and logging of the leak-tightness of systems containing hydrofluorocarbons in accordance with Regulation EC842/2006, incl. test book	05	DL00-0034
Maintenance	One-off maintenance service in accordance with maintenance schedule. Visual inspection of mechanical and electrical components, testing of all key functions. Calibration of a test temperature specified by the user in center of usable space without certificate	05	DL20-0500
SILVER 3-year maintenance contract	Maintenance service as contractually agreed, visual inspection of mechanical and electrical components, check of control response, 20% discount on spare parts, testing of all key functions, calibration of one test temperature specified by the user in the center of the usable space, without certificate	05	DL20-0830
Calibration services			
Temperature and humidity calibration	Expansion –Temperature and humidity calibration with 1 measuring point in center of chamber with 1 specified pair of values, including certificate	03, 04, 05	DL30-0302
	Temperature and humidity calibration with 1 measuring point in center of chamber with 1 specified pair of values, including certificate	03, 04, 05	DL30-0301
Temperature and humidity measurement according to DIN12880	Temperature measurement in accordance with DIN 12880 with 27 temperature measuring points and 1 humidity measuring point in center of chamber with a pair of values specified by the user, including certificate	03, 04, 05	DL30-0427
Temperature and humidity measurement, 18-1 measuring points	Temperature measurement with 18 temperature measuring points and 1 humidity measuring point in center of chamber with a pair of values specified by the user, including certificate	03, 04, 05	DL30-0318
Temperature and humidity measurement, 27-1 measuring points	Temperature measurement with 27 temperature measuring points and 1 humidity measuring point in center of chamber with a pair of values specified by the user, including certificate	03, 04, 05	DL30-0327
Temperature and humidity measurement, 9-1 measuring points	Temperature measurement with 9 temperature measuring points and 1 humidity measuring point in center of chamber with a pair of values specified by the user, including certificate	03, 04, 05	DL30-0309
Validation services			
Execution of IQ/OQ	Execution of IQ/OQ in accordance with qualification folder	05	DL42-0300
Execution of IQ/OQ/PQ	Execution of IQ/OQ/PQ in accordance with qualification folder	05	DL44-0500
Warranty service			
1-year warranty extension	The warranty is extended by 1 year from the delivery date, wear parts are excluded	-	DL50-0040

\* Notes > See last page



#### NOTES

- Condensation may occur in the area around the access port. Access ports may be placed in custom locations for an additional charge. UL mark is not granted when this option is used. 01
- 02
- Sensor calibration is performed in an accredited calibration laboratory. 03
- Calibration is performed according to the BINDER factory standard. 04
- 05 Quoted prices do not include travel costs. Please refer to the chapter on BINDER Service for travel costs for your region. Quoted prices for services performed in Switzerland do not include a country-specific added fee (available on request).

#### **BINDER GmbH**

Tuttlingen, Germany TEL +49 7462 2005 0 FAX +49 7462 2005 100 info@binder-world.com www.binder-world.com

#### BINDER Asia Pacific (Hong Kong) Ltd.

Kowloon, Hong Kong, P.R. China TEL +852 39070500 FAX +852 39070507 asia@binder-world.com www.binder-world.com

#### **BINDER Environmental Testing**

Equipment (Shanghai) Co., Ltd. Shanghai, P.R. China TEL +86 21 685 808 25 FAX +86 21 685 808 29 china@binder-world.com www.binder-world.com

#### **Representative Office for CIS**

Moscow, Russia TEL +7 495 988 15 16 FAX +7 495 988 15 17 russia@binder-world.com www.binder-world.com

#### **BINDER Inc.**

Bohemia, NY, USA TEL +1 631 224 4340 FAX +1 631 224 4354 usa@binder-world.com www.binder-world.us